



KJM 9250

bfprf1bfprf2 experiments on the AVI-600 spectrometer

Version 7.3

Topspin 1.3 Windows XP AVI600



© Professor Emeritus Alistair Lawrence Wilkins,
University of Waikato, New Zealand.
January 2020

bfprf1bfprf2 Experiments on the AVI-600

1.0 Introduction

Bruker pp's with CW presaturation on F1 require that it is applied at O1 = the spectral window midpoint. In cases where target compound signals are in the range 0-6 ppm and the HOD or the other line to be presaturated occurs in the vicinity of 5 ppm, this requires that the presaturation experiment is set up with O1 = 5 ppm and SW = 10 ppm.

*aw coded bfprf1bfprf2 experiments can be used to apply dual CW presaturation at **CNST9 Hz on F1** and at **CNST21 Hz on F2**, independent of one's choice of SW and O1.*

All power levels, pulse types and pulse times used in aw coded bfprf1bfprf2 2D experiments are by default prosol Table linked and no FQLISTS are required.

Optimum outcomes are achieved in aw coded bfprf1bfprf2 experiments when the default prosol Table linked F1 and F2 presaturation powers (**PL9** and **PL21** respectively) are adjusted to be sufficient to suppress target signals while not suppressing correlations arising from other nearby signals.

Depending on the proximity of genuine signals this may require the **PL9** and **PL21** power levels to be reduced by 3-12 db from their default prosol Table values. This can be achieved by reducing (*ie* attenuating) the **PL9** and **PL21** power levels by adding 3-12 db to their default prosol Table power levels.

Prior to setting up a **bfprf1bfprf2** experiment determine the frequencies in **Hz** of the two signals to be presaturated in a standard ¹H spectrum and ascertain the desired 2D-spectral window (**SW**) in **ppm** and its intended **O1 value** in **Hz** (= spectral window midpoint)

2.0 bfprf1bfprf2 Experiments

The following **bfprf1bfprf2** parameter sets are available on the **AVI-600**.

- 2.1 **cosybfprf1bfprf2**
- 2.2 **toscybfprf1bfprf2**
- 2.3 **clean-tocsybfprf1bfprf2**
- 2.4 **dipsi2bfprf1bfprf2**
- 2.5 **noesybfprf1bfprf2**
- 2.6 **roesybfprf1bfprf2** (with CW spin lock)
- 2.7 **roesy2bfprf1bfprf2** (with pulsed spin lock)

2.1 Cosybfprf1bfprf2 with CW presaturation on F1 at cnst9 Hz and F2 at cnst21 Hz

Parameter set: **awcosybfprf1bfprf2 (+ getprosol)**

Pulse programme: **awcosygpqfbfprf1bfprf2**

Type **eda** (enter) and enter **SW(F2) in ppm**, note the spectral window in **Hz** that appears in the **SWH(F2)** box and copy and paste this value into the **SWH(F1)** box.

Check **SWH(F2) = SWH(F1)** in Hz including all dp's.

Enter **O1 = O2 = CNST1 in Hz** (*in this order!*) where **O1** = the spectral window midpoint. Enter **CNST9** and **CNST21 in Hz** of the signals to be presaturated on F1 and F2\.

TD(F2) = 2K, **TD(F1) = 128-256** (your choice).

P0 = θ degree excitation pulse time, typically use a 45° or 90° pulse.

NS = 2, 4, 8 (any number is OK), **DS** = 2, 4 or 8.

D1 = presaturation time = 1.5, 2.0 sec or other time of your choice.

Type **ased** (enter) and review parameters used in the job including **CNST1**, **CNST9** and **CNST21**. Check gradient and **PL9** and **PL21** presaturation power levels are OK and if required adjust them appropriately.


Set receiver gain using **RGA** (*important!*).

Process with: **SI(F2) = SI(F1) = 2K**

WDW(F1) = WDW(F2) = SINE

SSB(F2) = SSB(F1) = 0

xfb, abs1, abs2 and (optionally) **sym**

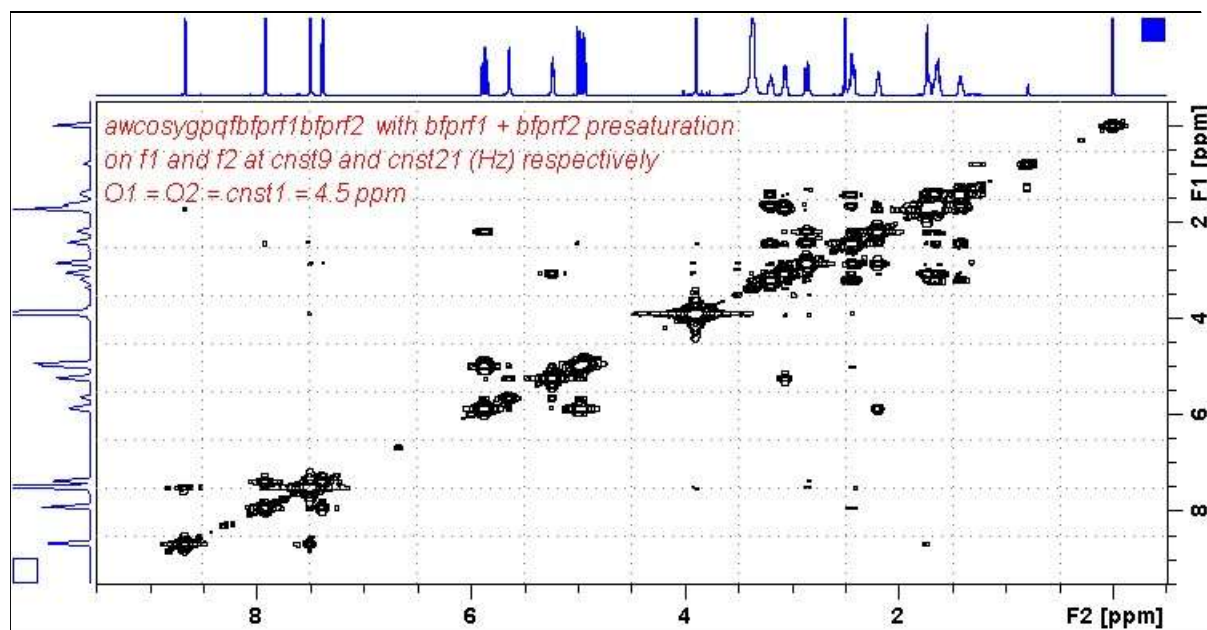

600 MHz **Cosybfprf1bfprf2** spectrum of quinine in D_6 -DMSO with the HOD and DMSO lines at 3.38 ppm and 2.5 ppm suppressed. The spectrum is centered at 4.5 ppm.

2.2 Tocsybfprf1bfprf2 with CW presaturation on F1 at cnst9 Hz and F2 at cnst21 Hz

Parameter set: **awtocsybfprf1bfprf2 (+ getprosol)**

Pulse programme: **awtocsybfprf1bfprf2**

Type **eda** (enter) and enter **SW(F2) in ppm**, note the spectral window in **Hz** that appears in the **SWH(F2)** box and copy and paste this value into the **SWH(F1)** box.

Check **SWH(F2) = SWH(F1)** in Hz including all dp's.

Enter **O1 = O2 = CNST1 in Hz (in this order!)** where **O1** = the spectral window midpoint.

Enter **CNST9** and **CNST21 in Hz** of the signals to be presaturated on F1 and F2.

TD(F2) = 2K, **TD(F1) = 128-256** (your choice).

NS = 4, 8, 16 (multiple of 4 or 8 recommended), **DS = 4 or 8**.

D1 = presaturation time = 1.5, 2.0 sec or other time of your choice.

D9 = correlation time = **80 msec** or other time of your choice (6-240 msec).

Type **ased** (enter) and review parameters used in the job including **CNST1**, **CNST9** and **CNST21**. Check the **PL9** and **PL21** presaturation power levels are OK and if required adjust them appropriately.


Set **receiver gain** using **RGA (important!)**.

Process with: **SI(F2) = SI(F1) = 2K**

WDW(F1) = WDW(F2) = QSINE

SSB(F2) = SSB(F1) = 2

xfb, abs1, abs2 and optionally **syma**

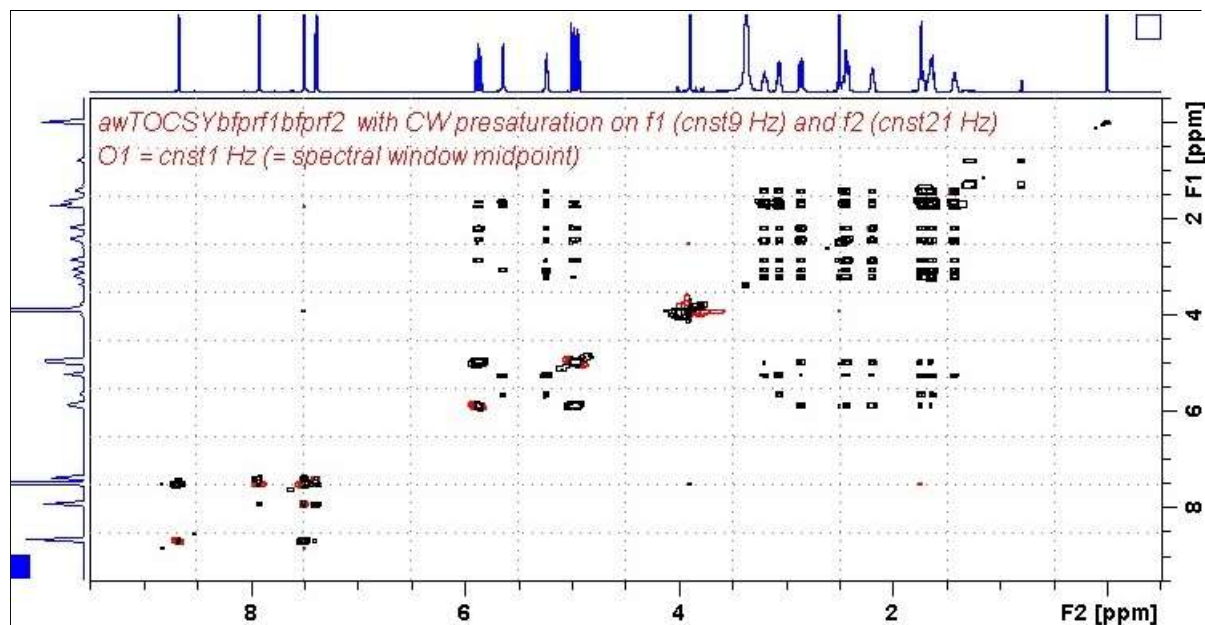

600 MHz **Tocsybfprf1bfprf2** spectrum of quinine in D₆-DMSO with the HOD and DMSO lines at 3.38 ppm and 2.5 ppm suppressed. The spectrum is centered at 4.5 ppm.

2.3 Clean-tocsybfprf1bfprf2 with CW presaturation on F1 at cnst9 Hz and F2 at cnst21 Hz

Parameter set: **awcleantocsybfprf1bfprf2 (+ getprosol)**

Pulse programme: **awclmlevbfprf1bfprf2**

Type **eda** (enter) and enter **SW(F2) in ppm**, note the spectral window in **Hz** that appears in the **SWH(F2)** box and copy and paste this value into the **SWH(F1)** box. Check **SWH(F2) = SWH(F1)** in Hz including all dp's.

Enter **O1 = O2 = CNST1 in Hz (in this order!)** where **O1** = the spectral window midpoint. Enter **CNST9** and **CNST21 in Hz** of the signals to be presaturated on F1 and F2.

TD(F2) = 2K, TD(F1) = 128-256 (your choice).

NS = 4, 8, 16 (multiples of 4 or 8 recommended), **DS = 4 or 8**.

D1 = presaturation time = 1.5, 2.0 sec or other time of your choice.

D9 = correlation time = 80 msec or other time of your choice (6-240 msec).

Type **ased** (enter) and review parameters used in the job including **CNST1, CNST9** and **CNST21**. Check the **PL9** and **PL21** presaturation power levels are OK and if required adjust them appropriately.


Set **receiver gain** using **RGA (important!)**.

Process with: **SI(F2) = SI(F1) = 2K**

WDW(F1) = WDW(F2) = QSINE

SSB(F2) = SSB(F1) = 2

xfb, abs1, abs2 and optionally **syma**

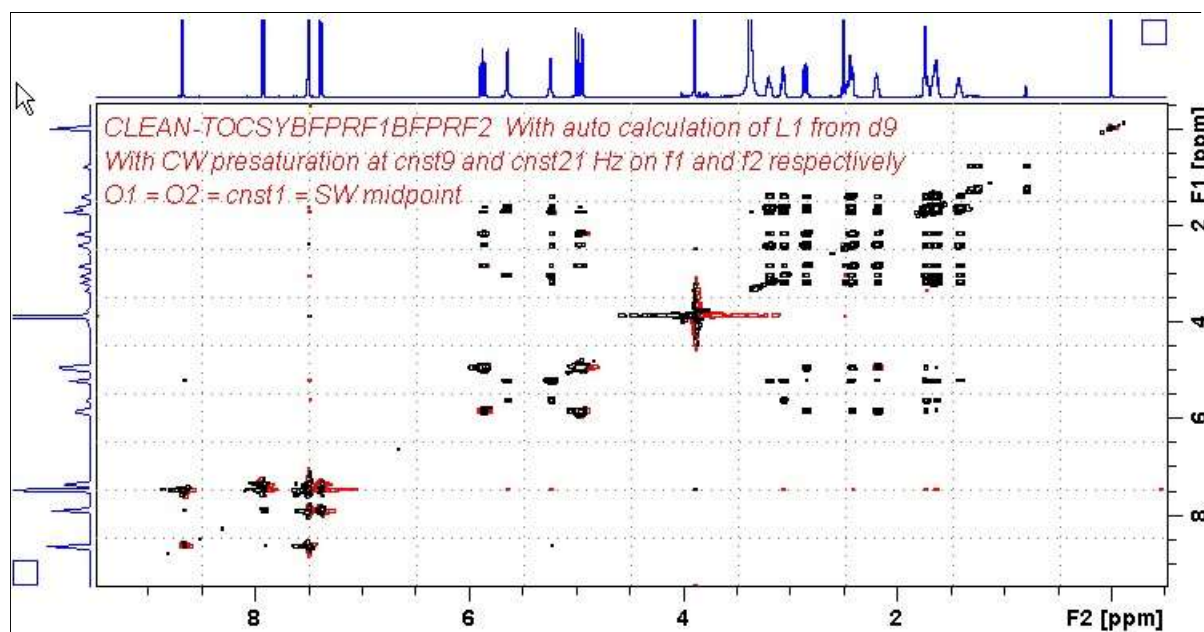

600 MHz **Clean-tocsybfprf1bfprf2** spectrum of quinine in D_6 -DMSO with the HOD and DMSO lines at 3.38 ppm and 2.5 ppm suppressed. The spectrum is centered at 4.5 ppm.

2.4 Dipsi2bfprf1bfprf2 with CW presaturation on F1 at cnst9 Hz and F2 at cnst21 Hz

Parameter set: **awdipsi2bfprf1bfprf2 (+ getprosol)**

Pulse programme: **awdipsi2phbfprf1bfprf2**

Type **eda** (enter) and enter **SW(F2) in ppm**, note the spectral window in **Hz** that appears in the **SWH(F2)** box and copy and paste this value into the **SWH(F1)** box. Check **SWH(F2) = SWH(F1)** in Hz including all dp's.

Enter **O1 = O2 = CNST1 in Hz** (*in this order!*) where **O1** = the spectral window midpoint. Enter **CNST9** and **CNST21 in Hz** of the signals to be presaturated on F1 and F2.

TD(F2) = 2K, **TD(F1) = 128-256** (your choice).

NS = 4, 8, 16 (multiples of 4 or 8 recommended), **DS = 4 or 8**.


D1 = presaturation time = 1.5, 2.0 sec or other time of your choice.

D9 = correlation time = **80 msec** or other time of your choice (6-240 msec).

Type **ascd** (enter) and review parameters used in the job including **CNST1**, **CNST9** and **CNST21**. Check the **PL9** and **PL21** presaturation power levels are OK and if required adjust them appropriately.

Set **receiver gain** using **RGA** (*important!*).

Process with: **SI(F2) = SI(F1) = 2K**
WDW(F1) = WDW(F2) = QSINE
SSB(F2) = SSB(F1) = 2
xfb, abs1, abs2 and optionally **syma**

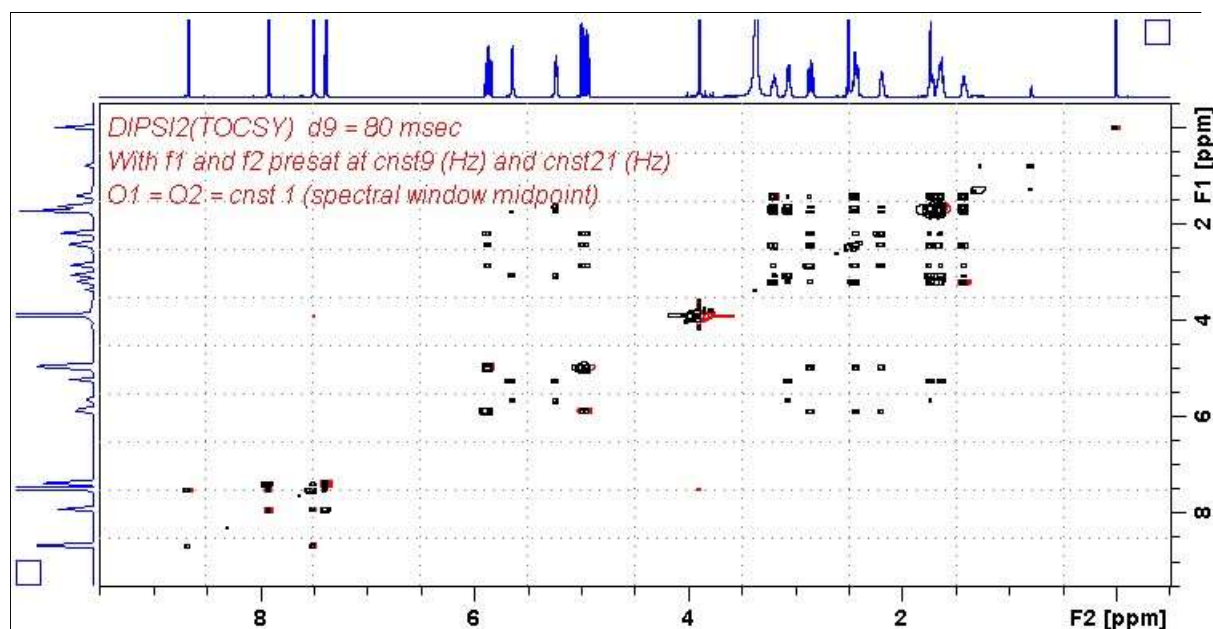

600 MHz **Dipsi2bfprf1bfprf2** spectrum of quinine in D_6 -DMSO with the HOD and DMSO lines at 3.38 ppm and 2.5 ppm suppressed. The spectrum is centered at 4.5 ppm.

2.5 Noesybfprf1bfprf2 with CW presaturation on F1 at cnst9 Hz and F2 at cnst21 Hz

Parameter set: **awnoesybfprf1bfprf2 (+ getprosol)**

Pulse programme: **awnoesygpphbfprf1bfprf2**

Type **eda** (enter) and enter **SW(F2) in ppm**, note the spectral window in **Hz** that appears in the **SWH(F2)** box and copy and paste this value into the **SWH(F1)** box.

Check **SWH(F2) = SWH(F1)** in Hz including all dp's.

Enter **O1 = O2 = CNST1 in Hz** (*in this order!*) where **O1** = the spectral window midpoint.

Enter **CNST9** and **CNST21 in Hz** of the signals to be presaturated on F1 and F2.

TD(F2) = 2K, **TD(F1) = 128-256** (your choice).

NS = 4, 8, 16 (multiples of 4 or 8 recommended), **DS = 4 or 8**.

D1 = presaturation time = 1.5, 2.0 sec or other time of your choice.

D8 = NOE mixing time = **0.5 sec** or other time of your choice.

Type **ased** (enter) and review parameters used in the job including **CNST1**, **CNST9** and **CNST21**. Check the **PL9** and **PL21** presaturation power levels are OK and if required adjust them appropriately.


Set receiver gain using **RGA** (*important!*).

Process with: **SI(F2) = SI(F1) = 2K**

WDW(F1) = WDW(F2) = QSINE

SSB(F2) = SSB(F1) = 2

xfb, abs1, abs2 and optionally **syma**

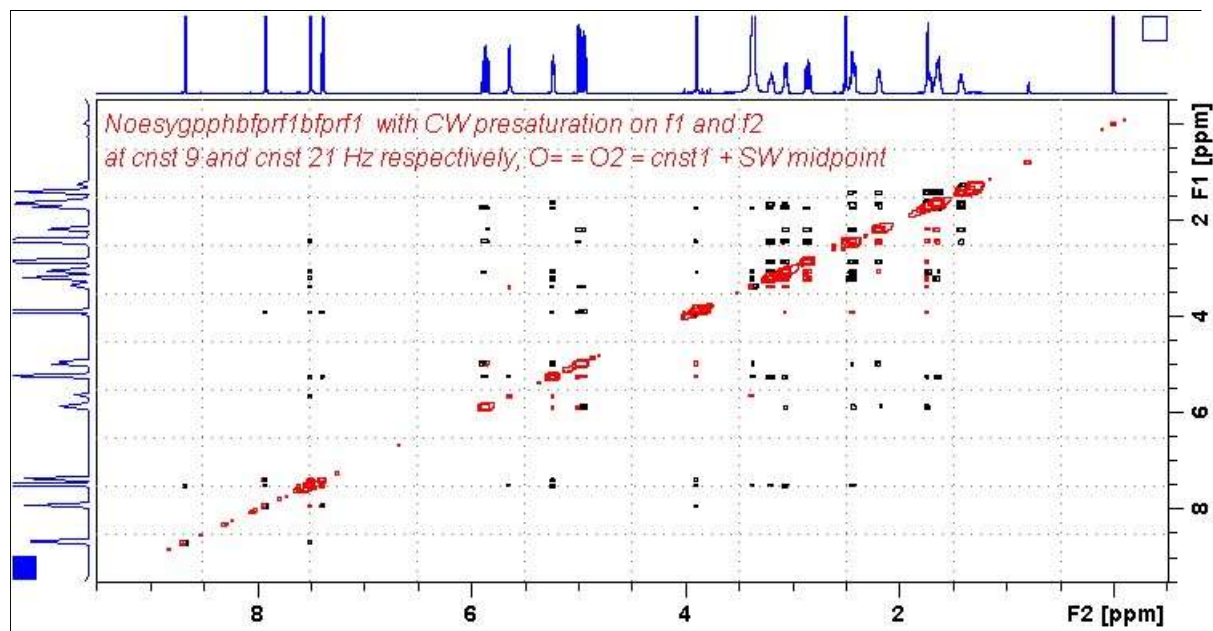

600 MHz **Noesybfprf1bfprf2** spectrum of quinine in D_6 -DMSO with the HOD and DMSO lines at 3.38 ppm and 2.5 ppm suppressed. The spectrum is centered at 4.5 ppm.

2.6 Roesybfprf1bfprf2 with CW presaturation on F1 at cnst9 Hz and F2 at cnst21 Hz

Parameter set: **awroesybfprf1bfprf2 (+ getprosol)**

Pulse programme: **awroesyphbfprf1bfprf2** (with CW spinlock)

Type **eda** (enter) and enter **SW(F2) in ppm**, note the spectral window in **Hz** that appears in the **SWH(F2)** box and copy and paste this value into the **SWH(F1)** box. Check **SWH(F2) = SWH(F1)** in Hz including all dp's.

Enter **O1 = O2 = CNST1 in Hz** (*in this order!*) where **O1** = the spectral window midpoint. Enter **CNST9** and **CNST21 in Hz** of the signals to be presaturated on F1 and F2.

TD(F2) = 2K, **TD(F1) = 128-256** (your choice).

NS = 4, 8, 16 (multiples of 4 or 8 recommended), **DS = 4 or 8**.


D1 = presaturation time = 1.5, 2.0 sec or other time of your choice.

P15 = spinlock time = **200000 usec** or **250000 usec** = 250 or 200 msec.

Type **ased** (enter) and review parameters used in the job including **CNST1**, **CNST9** and **CNST21**. Check the **PL9** and **PL21** presaturation power levels are OK and if required adjust them appropriately.

Set **receiver gain** using **RGA** (*important!*).

Process with: **SI(F2) = SI(F1) = 2K**
WDW(F1) = WDW(F2) = QSINE
SSB(F2) = SSB(F1) = 2
xfb, abs1, abs2 and optionally **syma**

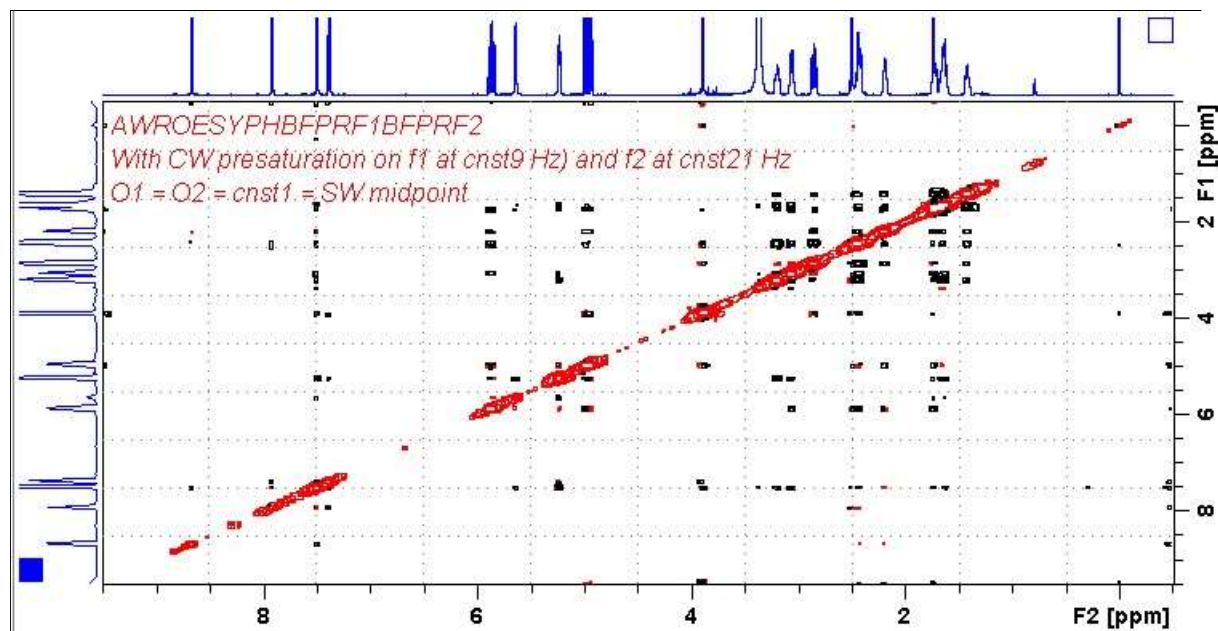

600 MHz **Roesybfprf1bfprf2** spectrum of quinine in D_6 -DMSO with the HOD and DMSO lines at 3.38 ppm and 2.5 ppm suppressed. The spectrum is centered at 4.5 ppm.

2.7 Roesy2bfprf1bfprf2 with CW presaturation on F1 at cnst9 Hz and F2 at cnst21 Hz

Parameter set: **awroesy2bfprf1bfprf2 (+ getprosol)**

Pulse programme: **awroesy2bfprf1bfprf2**(with pulsed spinlock)

Type **eda** (enter) and enter **SW(F2) in ppm**, note the spectral window in **Hz** that appears in the **SWH(F2)** box and copy and paste this value into the **SWH(F1)** box. Check **SWH(F2) = SWH(F1)** in Hz including all dp's.

Enter **O1 = O2 = CNST1 in Hz** (*in this order!*) where **O1** = the spectral window midpoint. Enter **CNST9** and **CNST21 in Hz** of the signals to be presaturated on F1 and F2.

TD(F2) = 2K, **TD(F1) = 128-256** (your choice).

NS = 4, 8, 16 (multiples of 4 or 8 recommended), **DS = 4 or 8**.

D1 = presaturation time = 1.5, 2.0 sec or other time of your choice.

P15 = spinlock time = **200000 usec** or **250000 usec** = 200 or 250 msec.

Type **ascd** (enter) and review parameters used in the job including **CNST1**, **CNST9** and **CNST21**. Check the **PL9** and **PL21** presaturation power levels are OK and if required adjust them appropriately.


Set **receiver gain** using **RGA** (*important!*).

Process with: **SI(F2) = SI(F1) = 2K**

WDW(F1) = WDW(F2) = QSINE

SSB(F2) = SSB(F1) = 2

xfb, abs1, abs2 and optionally **syma**

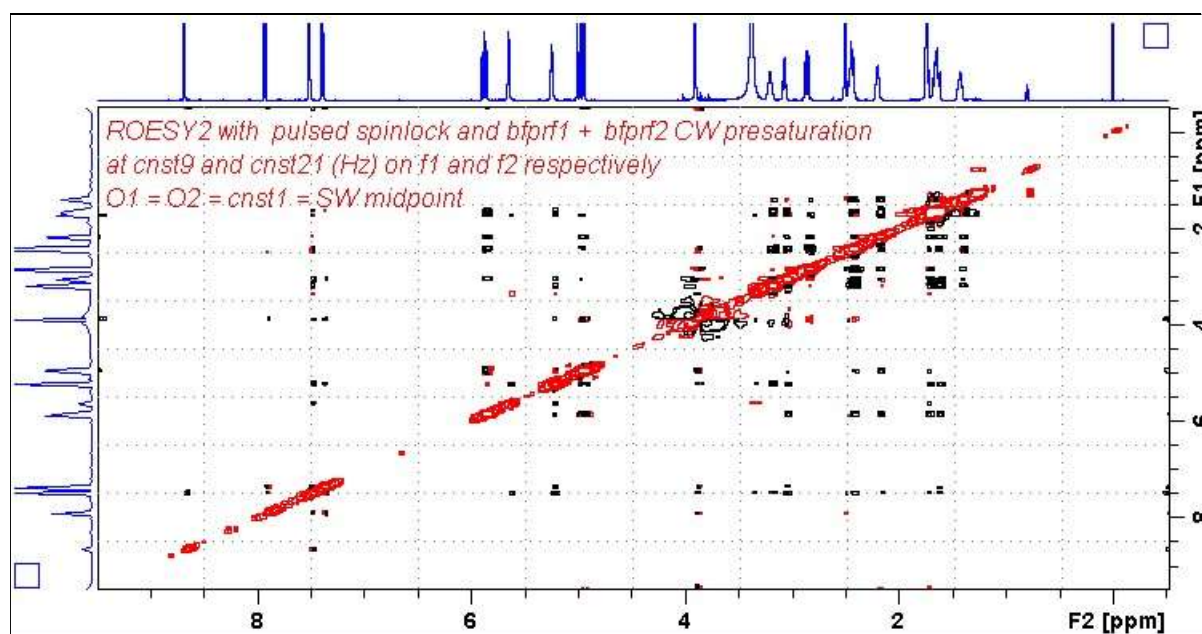

600 MHz **Roesy2bfprf1bfprf2** spectrum of quinine in D₆-DMSO with the HOD and DMSO lines at 3.38 ppm and 2.5 ppm suppressed. The spectrum is centered at 4.5 ppm.