




When you are connected to the UiO network click on Start in lower left corner and then on Computer. When connected to Eduroam and using vpn this might or might not work. Select Map network drive.

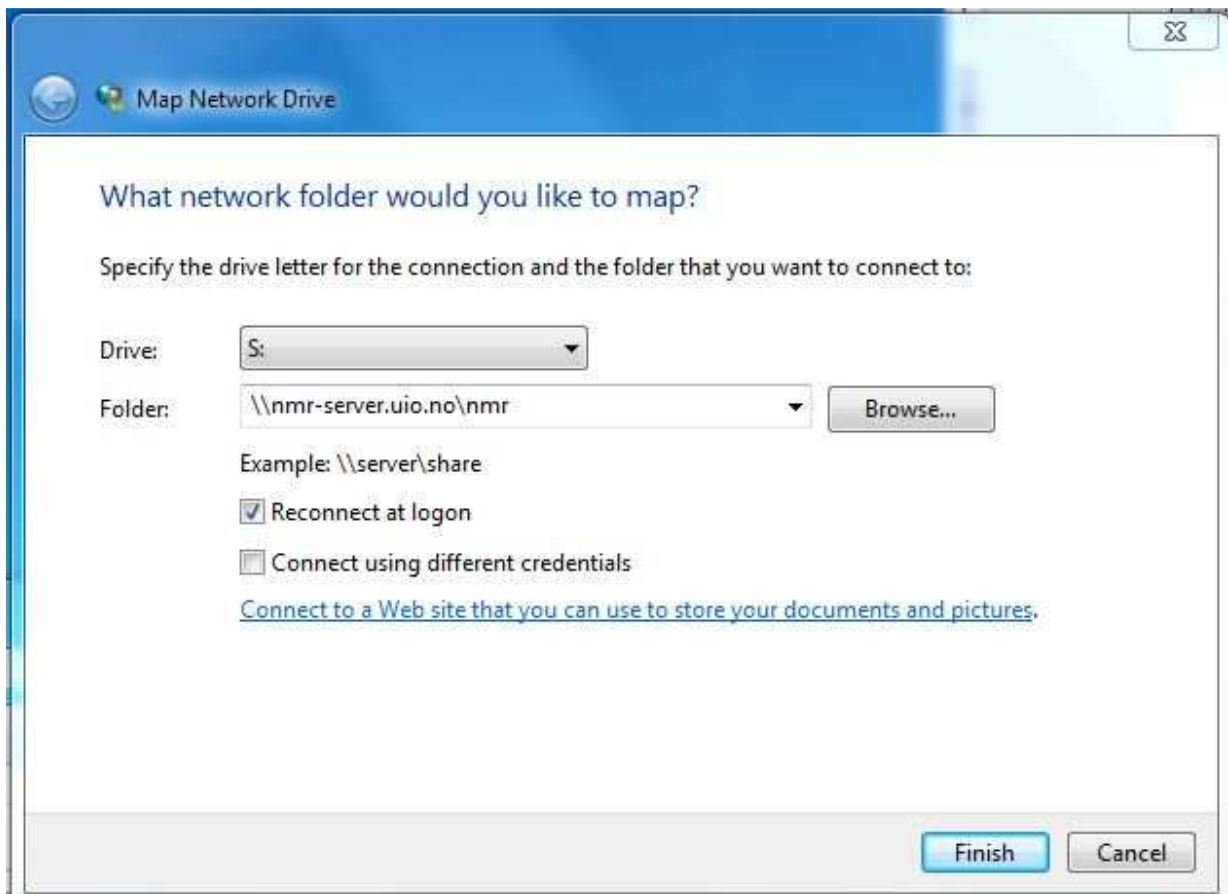

Type in as shown above. Click on Finish.

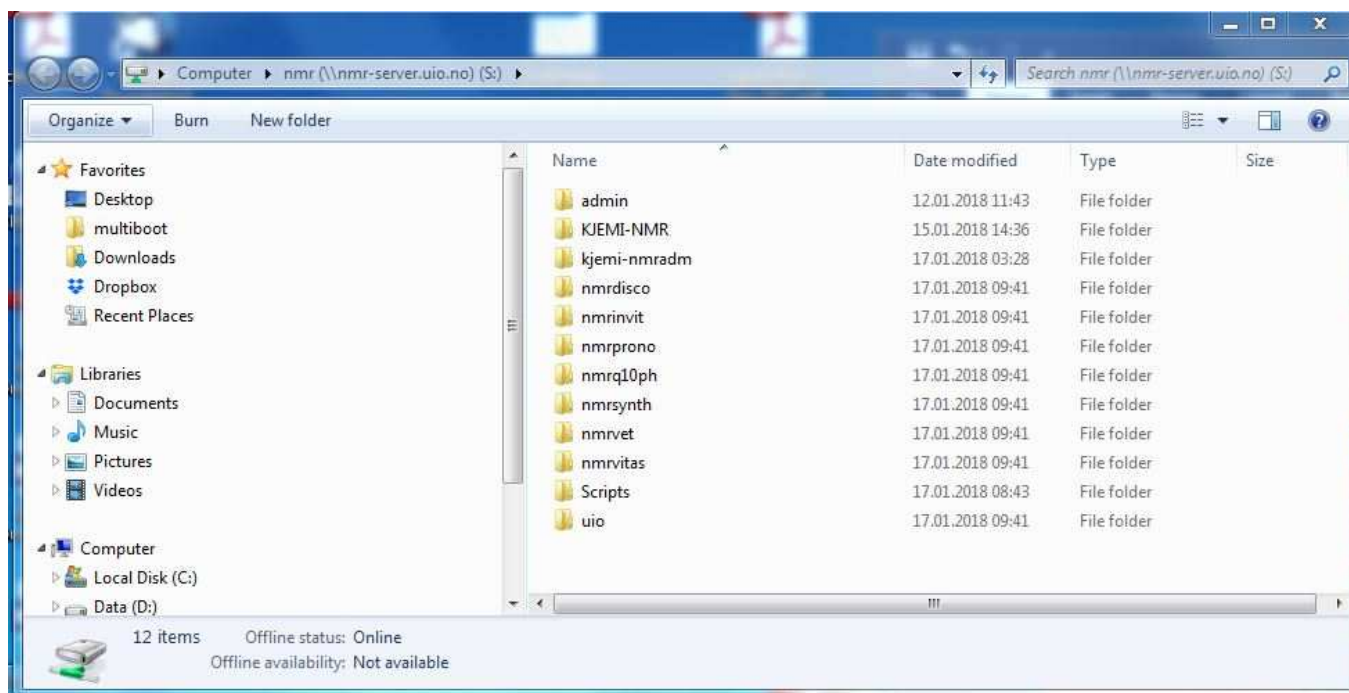

For most users locate the uio folder. Click on it.

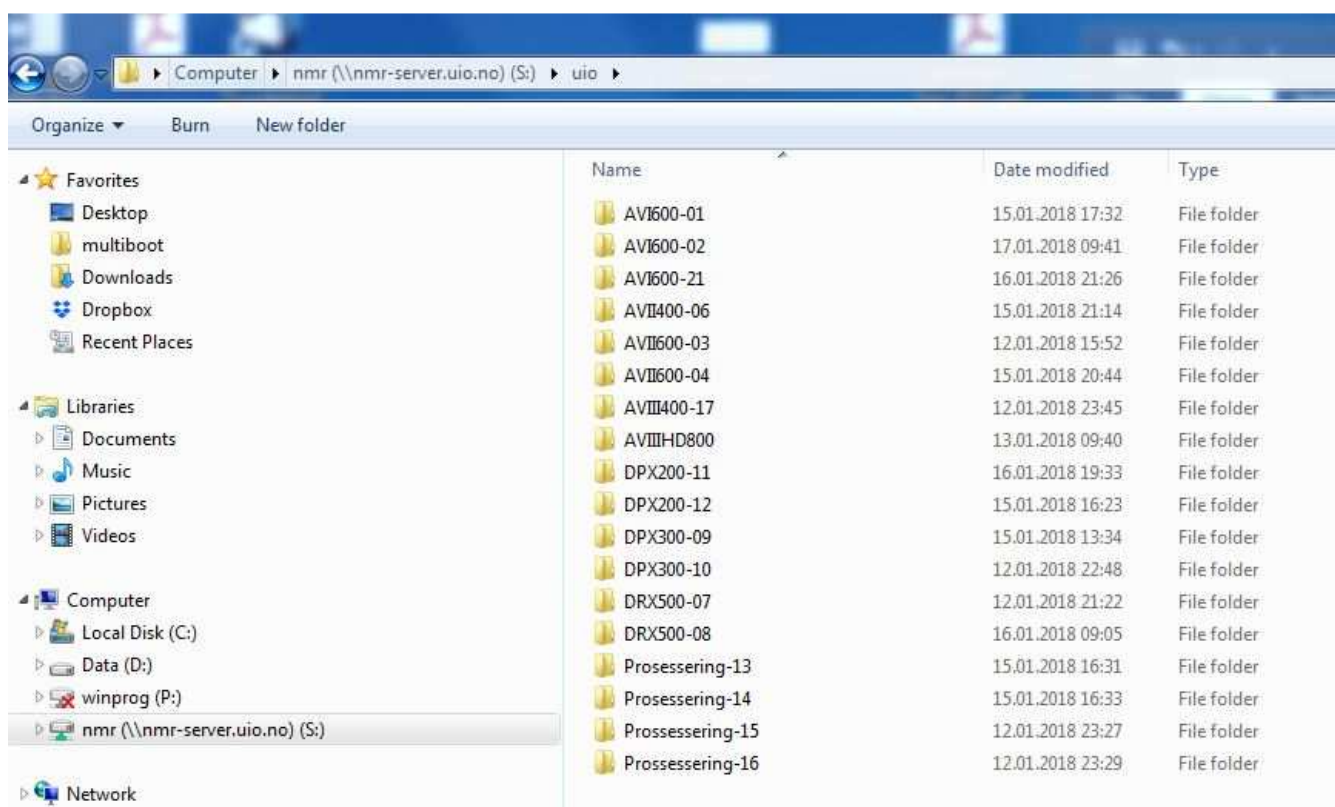

The nmr folders shows up. Some folders are missing here since the copying is not completed yet. Click on the relevant folder and all usernames will appear. Find yours and dig into it until you find your data. Drag and drop your nmr data to where you store them on your own PC or on the M drive. nmr-server.uio.no has 16 TB of storage space and data will not be deleted for many years to come. However, if the server burns out the data will be lost. Most likely it still reside on the nmr PCs where it originated. To be on the safe side you ought to make backup of new data every day.