



KJM 9250

**Proton $T_2$  Spectra on the DRX500 Spectrometer**

Version 5.0

Topspin 1.3 Windows XP DRX 500



© Professor Emeritus Alistair Lawrence Wilkins,  
University of Waikato, New Zealand.  
April 2018

## Proton T<sub>2</sub> Spectra on the DRX-500

### 1.0 Introduction

An **awprotont2** parameter set and linked **VCLIST** file from which a **VDLIST** file *must* be calculated prior to processing a T<sub>2</sub> data set have been set up on the DXR-500 spectrometer running under TS1.3.

The parameter sets have **D1 = 10 sec** and 8 x **VCLIST** values. The T<sub>2</sub> data set is acquired using the *smallest vclist value first. Prior to processing T<sub>2</sub> data sets variable constant (VC) values must be converted to time domain variable delay (VD) values using the **vctovd** command (see instructions 15 and 16.)*

**D1** and the longest **VD** time derived from the largest VC **value** should be 3-5 times the longest T<sub>2</sub> in the sample compound. T<sub>2</sub> is always shorter than T<sub>1</sub>.

$$VD_{\max} = [P2 + (2 * D20)] * VC_{\max}$$

T<sub>2</sub> data sets can be processed on the spectrometer terminal or off line using any version of Topspin provided the linked **VCLIST** file is either resident in the experiment's top level folder (= default setting: see below) or recreated or copied into the off line terminal's *C:\Bruker\TopspinX.X\exp\stan\nmr\lists\vc* folder.


If required the VCLIST file that is present in the experiment's top level folder can be opened with **WordPad** and the VC values in it can be viewed and used to recreate the original named **VCLIST** file from which the **VDLIST** file can be created using the **vctovd** command.

uxnmr.info	2/01/2017 10:50 p...	INFO File	4 KB
uxnmr.par	2/01/2017 10:50 p...	PAR File	24 KB
vclist	8/02/2018 10:30 p...	File	1 KB

NB: The experimental copy of the variable constant file is named as **vclist** irrespective of the name of the source vclist file.

## 2.0 Experiment Set Up

- 1) Create an experiment and read in the **awprotont2** parameter set (+ **getprosol**).  
Pulse programme = **cpmg**
- 2) Review default settings. These settings can be adjusted if required.  
**TD(F2) = 32K, TD(F1) = 8**  
**SW = 14 ppm, O1P = 6.3 ppm.**  
**D1 = 10 sec.**  
**D20 = 600 usec = 0.000600 sec**  
**NS = multiple of 2, 4 or 8, DS = 0, 2 or 4.**
- 3) Type **ased** (enter) and review other parameters used in the job including the linked **VCLIST** file = **AWPROTONT2-8**.
- 4) The **VCLIST** file should have the entries shown below, or other similar values. Do not alter the values in this file. An alternatively named **VCLIST** file should be created if different **VC** and **D1** values are required for a particular compound.

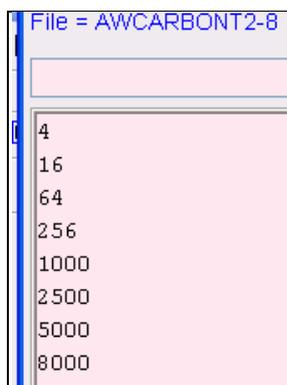


- 5) Set receiver gain using **RGA** (*important!*).
- 6) Type **edp** (enter) and check that **SI(F2) = 32K, SI(F1) = 8,**  
**WDW(F2) = EM, LB (F2) = 0.3-0.5 Hz** or other value of your choice.

SI	32768	8	Size of real spectrum
SF [MHz]	500.1300000	500.1300000	Spectrometer frequency
OFFSET [ppm]	13.30098	12.30805	Low field limit of spectrum
SR [Hz]	0	0	Spectrum reference frequency
HZpPT [Hz]	0.213709	751.201904	Spectral resolution
SPECTYP	UNDEFINED		Type of spectrum e.g. COSY, HI
Window function			
WDW	EM	no	Window functions for trf, xfb,...
LB [Hz]	0.30	0.30	Line broadening for em

- 7) Start the acquisition using the **ZG** command
- 8) When the experiment has run type **rser 1** (enter) to read in the first serial file which will appear in a TEMP screen display window.

9) Type **EFP** (return) to transform it and phase it as per a normal proton spectrum.

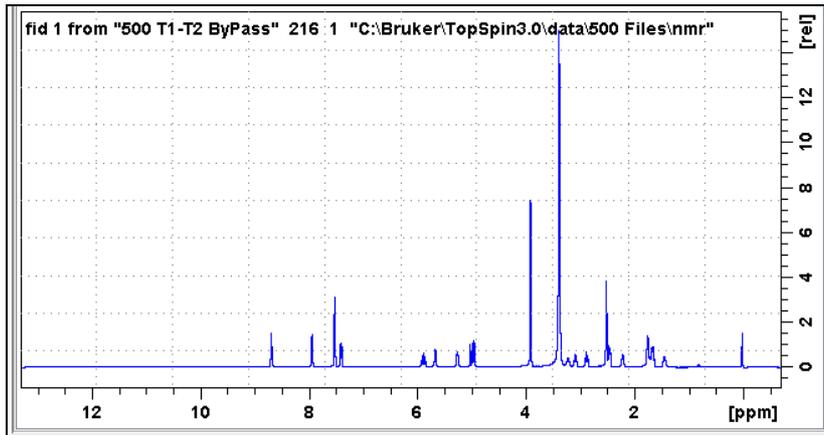

10) Type **edp** (enter) and note the phase constants for this spectrum.

PHC0 [degrees]	55.946
PHC1 [degrees]	-3.593
PH_mod	pk


11) Close the **TEMP** window and reload the T<sub>2</sub> data set file.

12) Type **edp** (enter) and enter the phase constants noted in step 10 above into the **F2 PHC0** and **F2 PHC1** cells and check **PH\_MOD** = pk. **F1** cell info is not used.

PHC0 [degrees]	55.946	0	0th order correction for
PHC1 [degrees]	-3.593	0	1st order correction for
PH_mod	pk	mc	Phasing modes for trf, x


13) Type **xf2** (return) to transform the 2D data set followed by **abs2** (return) to baseline smooth it.

14) Provided phase constants have been correctly set up the transformed data set plot should resemble that shown below. Signals intensities go from their maximum positive value to zero intensity.

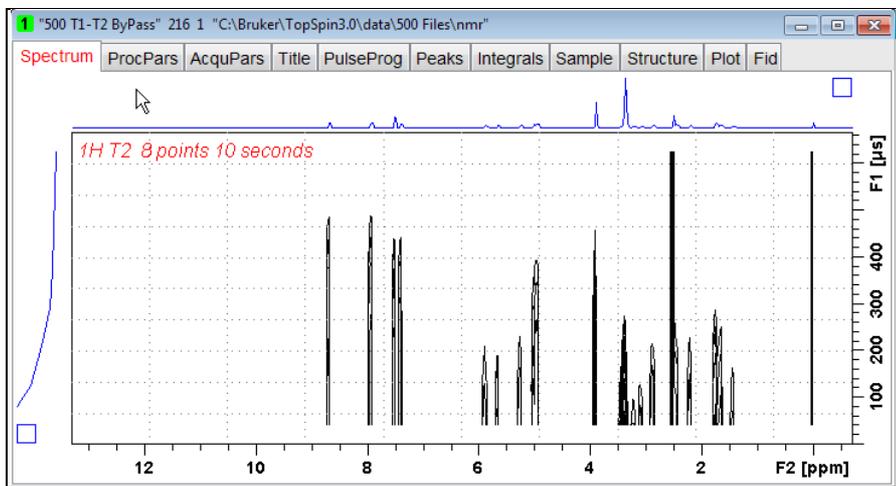


### 3.0 T<sub>2</sub> Data Set Processing

- 15) Prior to processing a T<sub>2</sub> data set it is essential that variable constant (VC) values are converted to time domain variable delay (VD) values.
- 16) Type **vctovd** (enter) and click OK on the display panels that appear.

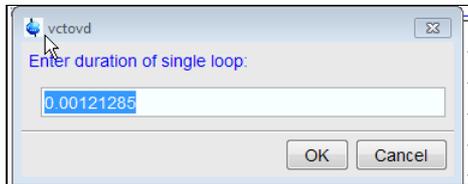


See the Appendix for the calculation of this value.

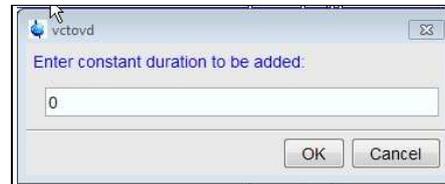


Not used.


- 17) Open Topspin's **Analysis** menu and click its **T1/T2 Relaxation** tab. If other Bruker processing software has been installed on the spectrometer terminal, as may be the case on the AVII-600, select the **Analysis** menu's **Topspin T1/T2 module** tab and open its **T1/T2 Relaxation** sub-menu tab.

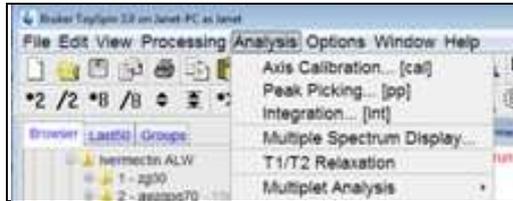


This will open up the screen display shown below.

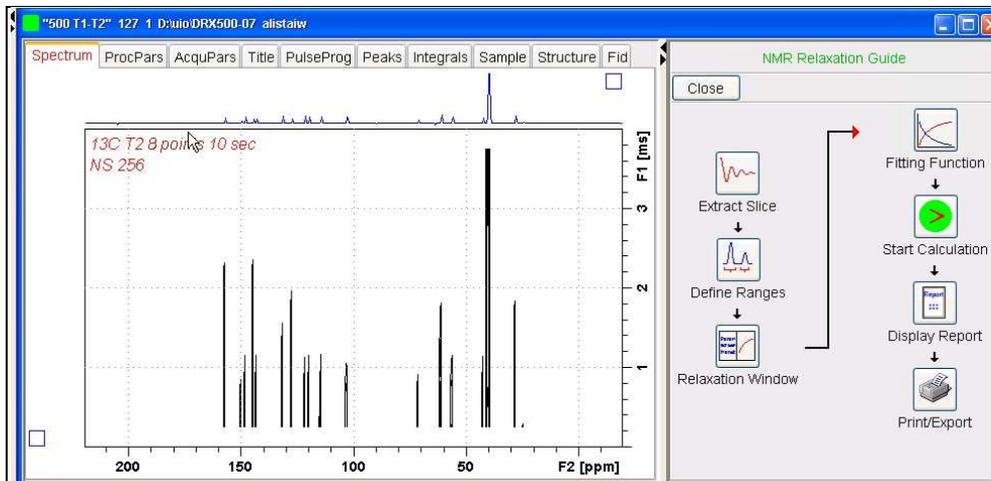


- 18) Click the **Extract Slice** button and then the click **Spectrum** button in the panel that appears.

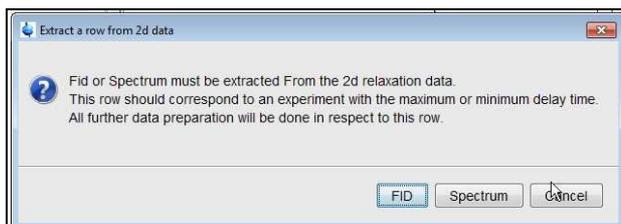


- 19) Enter **1** in the **Slice Number** cell and click the **OK** button. This will display the transformed spectrum ex the first (shortest) **VCLIST** → **VDLIST** value = the one that was phased via the **rser 1** routine in steps 8 and 9.

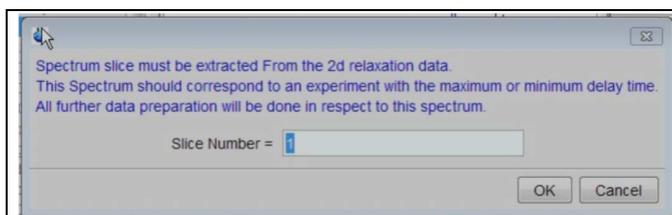

- 20) Expand the spectrum that appears in the usual way and integrate selected peaks in it. Integrals should start and terminate as close as possible to the edge of a peak.
- 21) Click the "**Save Region As**" icon button (= 7th from the left in the upper menu bar: *the one with the floppy disk icon*) and then its **Export Regions to Relaxation Module and .ret.** tab.

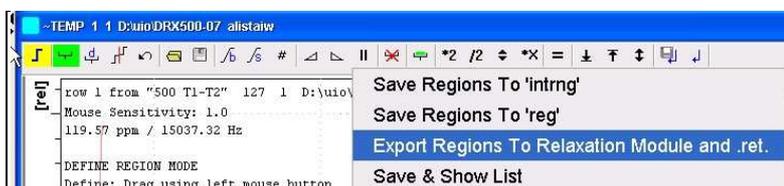


- 22) Click the **Relaxation Window** button and select **Intensity** or **Area** in the plot window that appears. **Area** is usually the better choice.

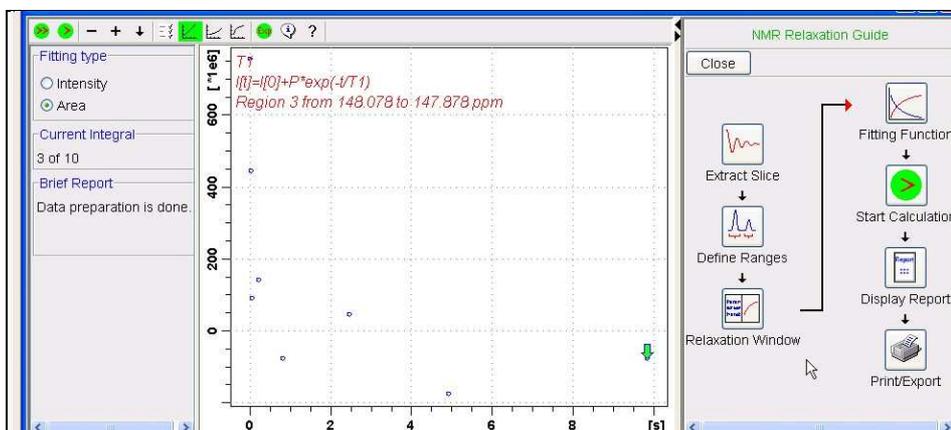


*The appearance of T1 information in this screen is a long standing Topspin bug.*

- 23) Click the **Fitting Function** button and note the comments about **intensity** or **area** options in the screen display that appears and close it.

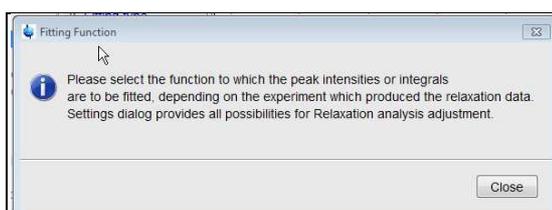


- 24) When the Fitting Function message screen is closed a panel with relaxation parameters (split into two screen captures below) will appear.

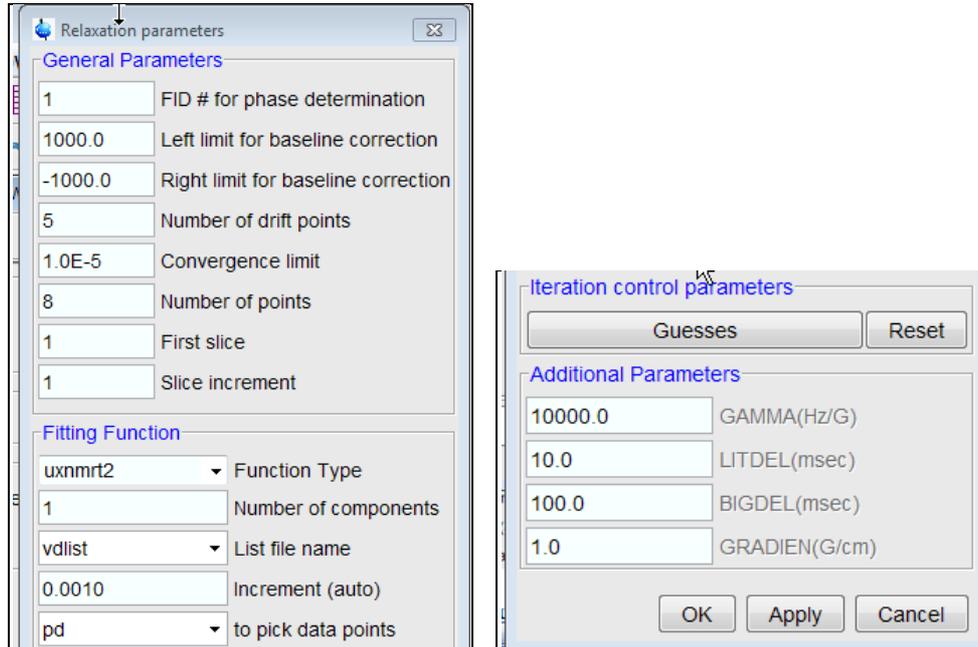

- 25) Check **Function Type = uxnmrt2** (*important!*) and **List file name = vclist**. Other cells/values can be left as they are (= default settings). Click **Apply** and **OK** and close the window.

- 26) Click the **double red arrow** in the menu bar at the top left of the plot window and **NOT** the single red Start Calculation arrow button in the NMR Relaxation Guide menu below the Fitting Function button.

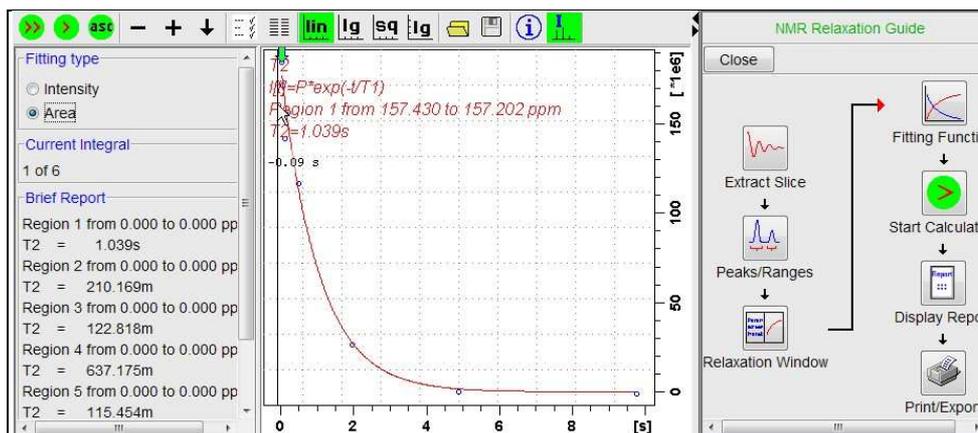

*The appearance of the  $T_1$  equation in this screen is a long standing Topspin error.*

- 27) The + and - buttons in the upper menu bar can be used move through the series of  $T_2$  plots.
- 28) Fitting type (**Intensity** or **Area**) can be changed in the plot display. If this is done clicking the **double red arrow** will recalculate the  $T_2$  results and update their plots.

29) The NMR Relaxation Guide has buttons which can be used to display and/or print T<sub>2</sub> results. A sample report for one signal is shown below. Zero value points (= completely T<sub>2</sub> relaxed) will be eliminated.

1 Dataset :  
 2 C:\Bruker\TopSpin3.0\data\500 Files\nmr\500 T1-T2 ByPass\216/pdata/1  
 3 AREA fit :  
 4  $I[t] = P \cdot \exp(-t/T_2)$   
 5  
 6 8 points for Integral 1, Integral Region from 0.000 to 0.000 ppm  
 7 Results Comp. 1  
 8  
 9 P = 9.147e-001  
 10 T2 = 1.994s  
 11 SD = 4.210e-002  
 12  
 13 tau ppm integral intensity  
 14  
 15 4.851m 8.680 1.1457e+009 2.7347e+007  
 16 19.406m 8.680 1.0785e+009 2.6697e+007  
 17 77.622m 8.680 9.7667e+008 2.3493e+007  
 18 310.490m 8.680 8.4007e+008 1.9537e+007  
 19 1.213s 8.680 5.5076e+008 1.2981e+007  
 20 3.639s 8.679 2.1239e+008 5.0751e+006  
 21 7.277s 8.679 4.7105e+007 1.1544e+006  
 22 14.554s 8.680 7.0986e+006 2.5912e+005

