

KJM 5280




MestreNova

Program installation and Introductory use


Version 1.1

© F. Rise May 2018

Please locate mestrelab.com/download/mnova/

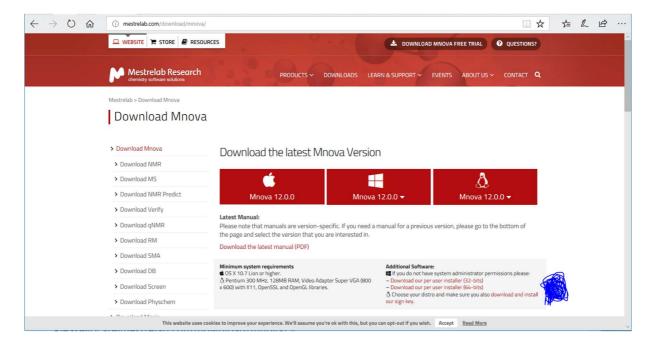

You are using Macintosh? Windows? Alternatively, Linux? When using Windows please look to the left of the blue clutter in the picture above. Download your version. In the following, the "Download per user installer (64-bits)" was used.

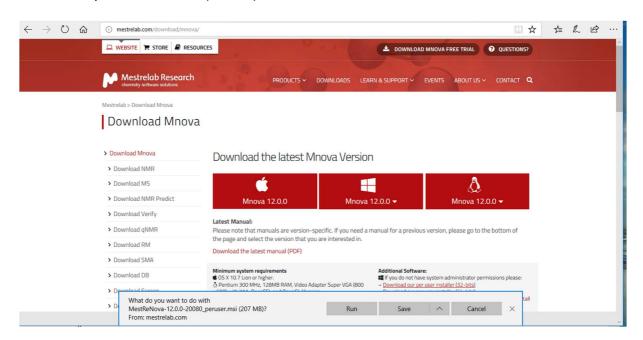


Save to desktop so you find what you downloaded.




Click on it.




Well? Next.

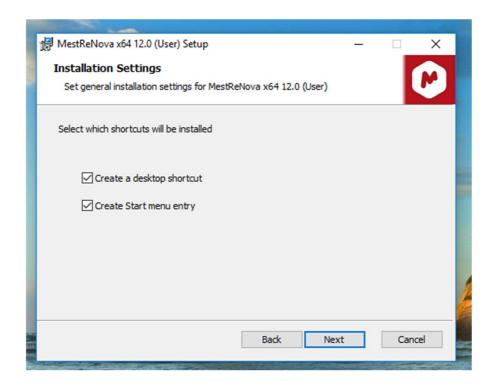


Next again.

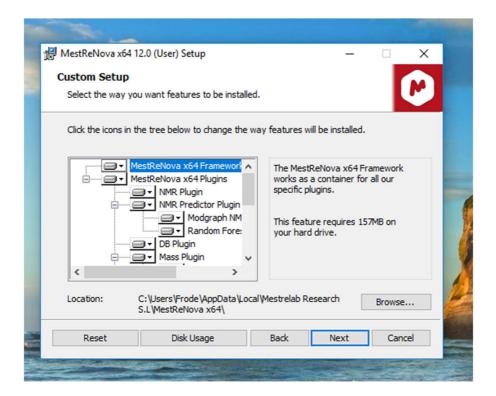


Whatever. No one understand this. Next.

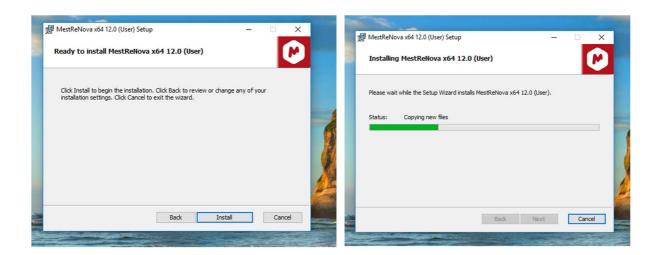

So many steps. Go on.

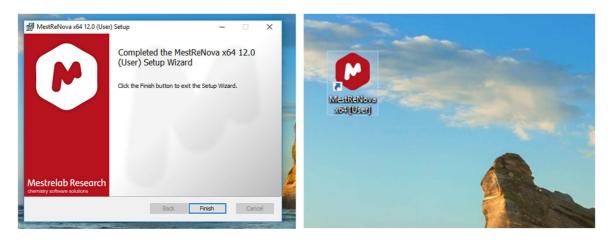

Finally, the program symbol/icon is on desktop. Click on it.

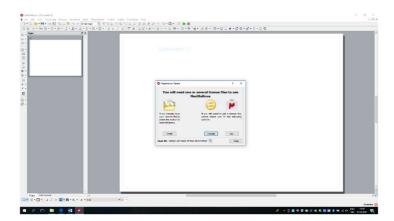


Oh – do I need a license? As a UiO user at either the Department of Chemistry or School of Pharmacy, you will obtain one, or rather three of them from either itsupport@kjemi.uio.no when

you send the physical address/MAC address of your PC/Mac to itsupport@kjemi.uio.no. Yes, your PC does indeed have a MAC address.


Go to Start in lower left corner of your Windows 7 PC. Type cmd in the command line.

In you type ipconfig -all


The physical address/MAC address in this instance is 56-A9-18-9D-AE-34 and not the IPv4 number.

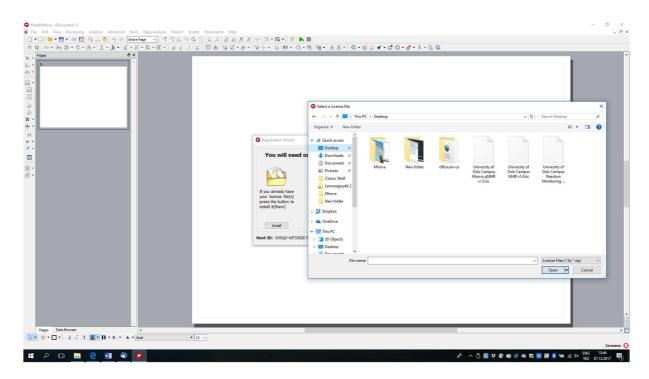


Locate your licenses. Click on install and be a little intelligent.

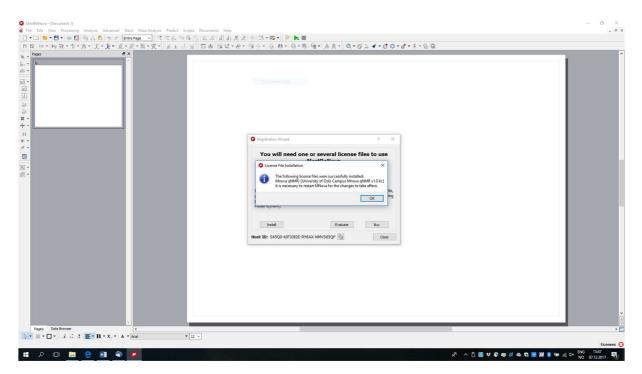

Install all three of them, one by one.

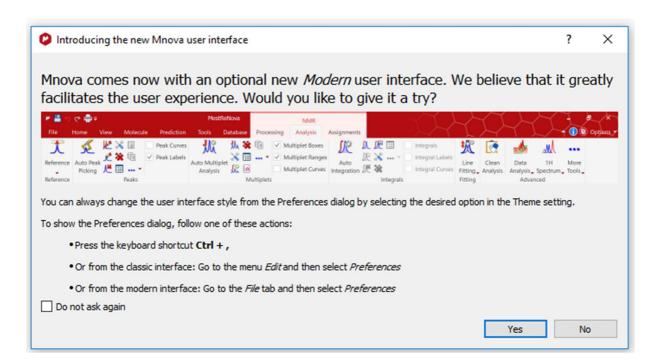


I really do not know what to suggest here. For the time being, the old user interface is used in this document. Restart MestreNova.

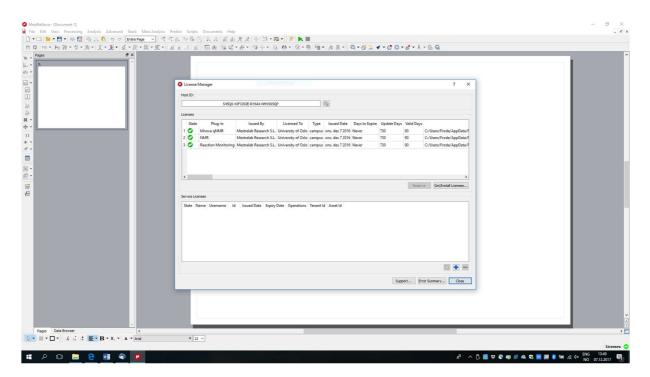


Help and License manager brings up this window. Note "Days to Expire" = Never. This is just as good as when you are collecting frequent flyer miles with Delta. SAS Euro bonus and KLM's customer program always let the miles expire for the author before they can be used for

anything. Fly Delta © and order your tickets for KLM destinations at https://www.delta.com/ as well.

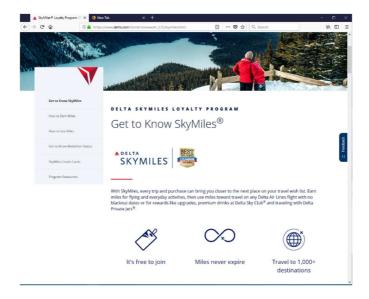


You can update MestreNova until 2019-12-07. The program license will never expire. When 150 users have installed MestreNova, the next user installing the program will not be able to use the program before one of 150 have forgotten to use her license for 90 days.

Before trying to explain intricacies of the program, the reader is advised to watch a bunch of YouTube videos showing how to use the program. Simply search for MestreNova and video. At the time, this document was written the following videos were instructive. Stop the videos often so you can repeat all the steps with your own nmr-data.

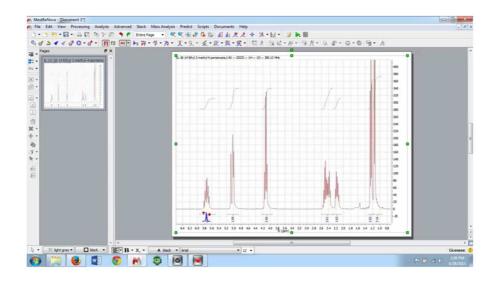


https://www.youtube.com/watch?v=DX02Nubdm44

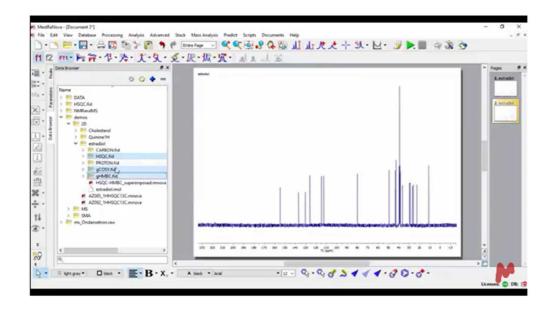

http://resources.mestrelab.com/mnova-2d-nmr-processing-a-quick-introduction/



https://www.youtube.com/watch?v=NqELnbaNI7s

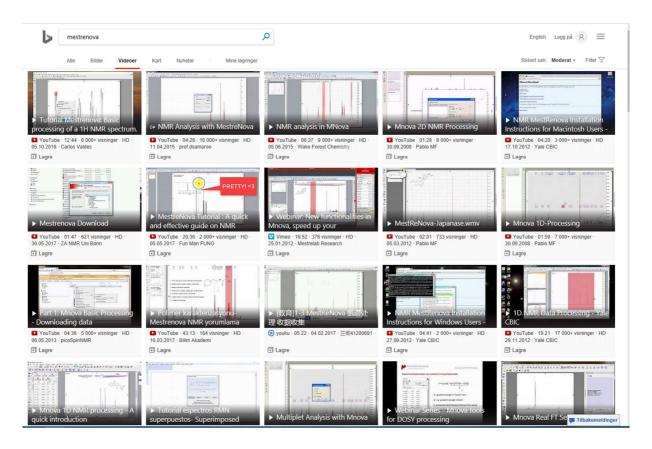

Searching the Internet brings up many MestreNova videos.